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


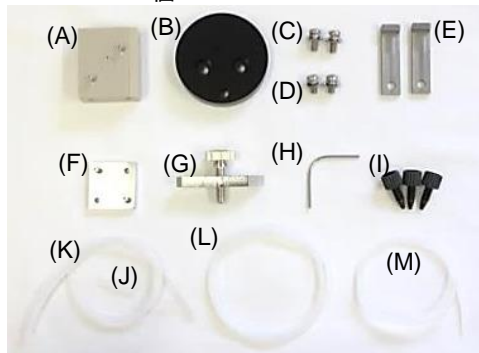
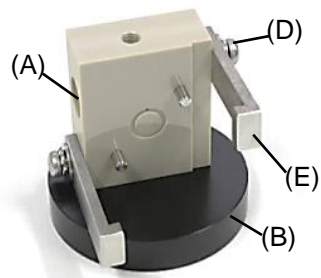
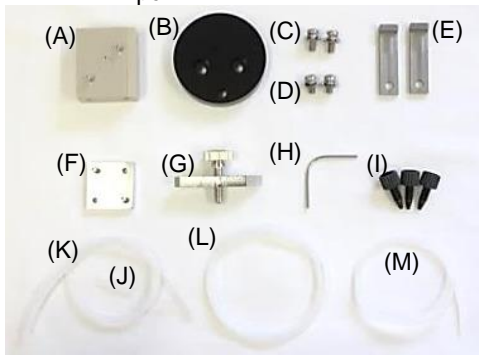
•Assembly

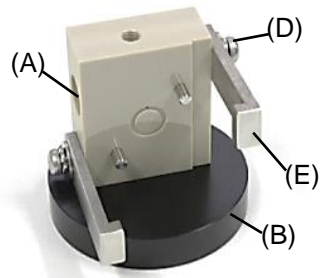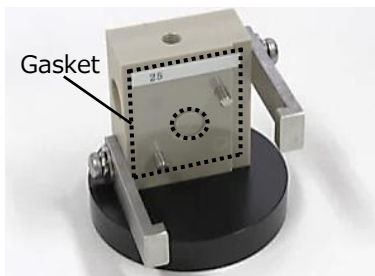
1. The radial flow cell block(A) is attached to the flow cell base block(B) with screws(C) from the bottom of the base, and flow cell arms(E) are connected to the electrode block with screws(D). Please use the Phillips head screwdriver for the assembly.

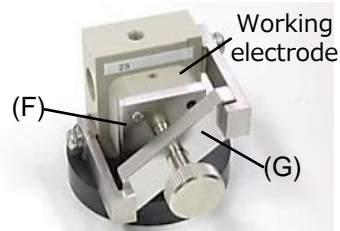
2. A gasket(option) is put on the electrode block as shown the figure to the left.

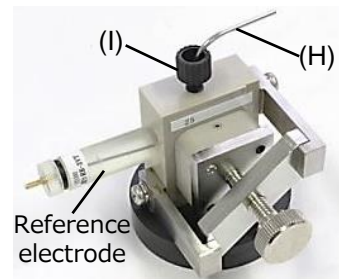3. The electrode surface of a working electrode(option) is put on the hole of the gasket,, and the stainless plate for clamp(F) is put and fixed as the left figure in the clamp(G).

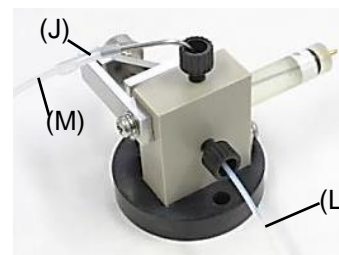4. A reference electrode(option), a fitting PEEK(I), and the stainless pipe 5cm(H) are attached as the left.

5. The Teflon tube(L) is connected with a fitting. Put the silicone tube(J) on the tip of the stainless pipe and connect the Teflon tube (M).

----- < Option > -----

- * Silicone tube(K) is a spare part of (J). You can also use it when you want to extend the outlet line, and so.
- * Don't tighten the clamp and screws too much when you set electrodes, fittings, and arms.
- * Put out the inlet tube from the tip of the fitting about 2mm when you connect them.

- Working Electrode
 - 012124 Glassy carbon electrode (Single 3mm)
 - 001016 Glassy carbon electrode (Single 6mm)
 - 000999 PFCE Carbon electrode (Single 3mm)
 - 011155 Gold electrode (Single 3mm)
 - 009908 Platinum electrode (Single 3mm)
 - 010251 Carbon paste electrode (Single 3mm)
- Reference Electrode
 - 013488 RE-3VT Reference electrode screw type (Ag/AgCl)
 - 013489 RE-7VT Non Aqueous reference electrode screw type (Ag/Ag+)
- Gasket
 - 001146 TG-2MR Teflon Gasket(Radial Flow)/12um (4pcs)
 - 001147 TG-5MR Teflon Gasket(Radial Flow)/25um (4pcs)
 - 001148 TG-6MR Teflon Gasket(Radial Flow)/50um (4pcs)
 - 012802 TG-8MR Teflon Gasket(Radial Flow)/100um(4pcs)

Please check color pdf data and the latest information with our website.

BAS Inc.

<https://www.als-japan.com> email: sales@als-japan.com