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<https://www.bas.co.jp> E-mail: sales@bas.co.jp

Working electrode manual

Thank you for your purchase from BAS Inc.

Please read following precautions before using products. If electrodes are used in violation of following precautions, any damage of the electrodes is not in our guarantee. Please check each serial code of label on electrodes, and search their pre-shipment check data from our website, and test electrodes under same experimental conditions to confirm initial quality of the electrodes.

(<https://www.als-japan.com/dl/>)

CAUTION

1. Do not remove each label on products or electrodes. The label is required for any case of inquiry.
2. Avoid using long time or intensive ultrasonic cleaning. Otherwise electrodes may possibly be damaged.
3. Do not use or wash electrodes with highly concentrated strong acidic or basic solutions.
4. Electrodes should be protected from strong shock.
5. Use electrodes under room temperature and atmospheric pressure.
6. Electrode modification by chemical or physical process will change electrode surface state. In this case, the electrode surface regeneration is not warranted.
7. Please use 013223 PK-3 Electrode Polishing kit for electrode polishing. Check the correct polishing method in our website.
(<https://www.als-japan.com/1735.html>)
8. Do not soak electrodes in organic solvents for long time, to avoid generating cracks in it.
9. Use electrodes on your own responsibility if you use instruments of other companies.
10. It is recommended that rotation rates of electrodes for RRDE-3A are not higher than 3000 rpm.
11. Do not disassemble working electrodes other than DRE disk replaceable electrodes.

note:

Initial defect due to our manufacturing (or production) process can be exchanged with a new product. Otherwise, problems happened by your experimental process or other reasons are not in our warranty.

Any defects in appearance causing no influence to measurement are not in our warranty.

Please check latest information on our website.

(<https://www.als-japan.com/support-product-manual.html>)

BAS Inc.

<https://www.als-japan.com> email: sales@als-japan.com